

Western Blot Analysis: Detection of Proteins using Chemiluminescence

Section of Cancer Genomics, Genetics Branch, NCI
National Institutes of Health

Reagents

ECL Western Blotting Analysis System

Amersham Pharmacia, Cat. RPN2108

Glycine

BioRad, Cat. 161-0716

Hyperfilm ECL

Amersham Pharmacia, Cat. RPN2103H

Methanol

Non-fat Dry Milk (NFDM)

Carnation Brand, Grocery Store of Choice

PVDF Membrane

BioRad, Cat. 162-0181

NaCl

Mallinckrodt, Cat. 7581

Tris Base

BioRad, Cat. 161-0717

Tris, pH 7.5, 1 M

Quality Biological, Inc., Cat.351-006-100

Tween 20

Sigma, Cat. P1379

Preparation

10X Running Buffer

Tris Base	30.3 g	f.c. [250 mM, pH 8.3]
-----------	--------	-----------------------

Glycine	144.2 g	f.c. [1.92 M]
---------	---------	---------------

Milli-Q water	1.0 L	
---------------	-------	--

Store at 4°C

1X Running Buffer

10X Running Buffer	100 ml	
--------------------	--------	--

Milli-Q Water	890 ml	
---------------	--------	--

10% SDS	10 ml	f.c. [0.1%]
---------	-------	-------------

Transfer Buffer

Tris Base	3.03 g	f.c. [25 mM, pH 8.3]
Glycine	14.4 g	f.c. [192 mM]
Milli-Q water	1.1 L	

10X TTBS

1M Tris (pH 7.5)	500 ml	f.c. [0.5 M]
NaCl	87.66 g	f.c. [1.5 M]
Tween-20	10 ml	f.c. [1%]

*Bring volume to 1L with sterile distilled water

Store at 4°C

1X TTBS

10X TTBS	100 ml
Milli-Q Water	900 ml

Store at 4°C

Procedure

1. Run Western gels in pre-chilled 1X Running Buffer at 200V for 60 min in ice bucket.
2. Transfer to PVDF membrane
 - a) Pre-wet PVDF in 50% MeOH starting at one end and slowly proceeding across membrane.
 - b) Equilibrate in 1X Transfer Buffer ~ 5 min.
 - c) Set-up transfer apparatus.
 - d) Transfer in pre-chilled 1X Transfer Buffer at 100 V for 60 min with ice pack and apparatus in ice bucket during transfer.
3. Block membrane in 20 ml 5% NFD, 1X TTBS for 60 min on shaker.
4. Detect membrane in 2 ml (1° Ab diluted in 5% NFD, 1X TTBS) in sealed bag on rocker overnight.
5. Rinse membrane 3 x 5 min with 1X TTBS.
6. Detect membrane in 10 ml (2° HRP-conjugated Ab + 1xTTBS) for 60 min, shaking
7. Rinse membrane 1 x 15 min followed by 4 x 5 min with 1X TTBS.
8. Mix equal volume Detection solutions 1 & 2 (125 µl/cm² membrane).

9. Add detection mixture to membrane allowing surface tension to keep solution on membrane surface without spilling off sides.
10. Incubate for exactly 1 min at RT.
11. Drain off detection solution and wrap membrane in Saran Wrap, carefully removing air pockets.
12. Expose to Hyperfilm for desired length of time.